

SpreetaTM

Immobilization of Ligand: Method 1

Application Brief

Number 001

2002

Sensors & Controls

IMPORTANT NOTICE

Texas Instruments and its subsidiaries (TI) reserve the right to make changes to their products or to discontinue any product or service without notice, and advise customers to obtain the latest version of the relevant information to verify, before placing orders, that information being relied on is current and complete. All products are sold subject to the terms and conditions of sale supplied at the time of order acknowledgement, including those pertaining to warranty, patent infringement, and limitation of liability.

TO warrants performance of its semiconductor products to the specifications applicable at the time of sale in accordance with TI's standard warranty. Testing and other quality control techniques are utilized to the extent TO deems necessary to support its warranty. Specific testing of all parameters of each device is not necessarily performed, except those mandated by government requirements.

CERTAIN APPLICATIONS USING SEMICONDUCTOR PRODUCTS MAY INVOLVE POTENTIAL RISKS OF EDEATH, PERSONAL INJURY, OR SEVERE PROPERTY OR ENVIRONMENTAL DAMAGE ("CRITICAL APPLICATIONS"). TI SEMICONDUCTOR PRODUCTS ARE NOT DESIGNED, AUTHORIZED, OR WARRENTED TO BE SUITABLE FOR USE IN LIFE SUPPORT DEVICES OR SYSTEMS OR OTHER CRITICAL APPLICATIONS. INCLUSION OF TI PRODUCTS IN SUCH APPLICATIONS IS UNDERSTOOD TO BE FULLY AT THE CUSTOMERS RISK.

In order to minimize risks associated with the customers product design. TI does not warrant or represent that any license, either express or implied, is granted under any patent right, copyright, mask work right, or other intellectual property right of TI covering or relating to any combination, machine, or process in which such semiconductor products or services might be or are used. TI's publication of the information regarding any third party's products or services does not constitute TI's approval, warranty or endorsement thereof.

Copyright ©2002Texas Instruments Incorporated

Immobilization of Ligand: Method 1

ABSTRACT

In this method the ligand is linked to the gold surface via a cross-linked protein film using carbodiimide mediated amide bond formation.

Introduction

It is well documented that direct physisorption of protein to a gold surface induces denaturation of the adsorbed protein. Here we assemble a cross-linked protein film on the gold surface and then link the ligand of interest. Thus the ligand will not be in direct contact with the gold hence preserving its binding activity. Figure 1 is an illustration depicting the immobilization process while Figure 2 is an actual response curve for the immobilization of antibody using the method outlined below.

Materials & Methods

All reagents were obtained from Sigma-Aldrich unless otherwise stated. Note: Always use gloves when handling reagents and sensors. Ensure adequate safety precautions are taken when handling all reagents especially succinic anhydride, N-hydroxysuccinimide (NHS) or N-ethyl-N'(dimethylaminopropyl)-carbodiimide (EDC). Refer to relevant MSDS for guidelines.

Preparation of Cross-linked Carboxymethylated Protein Film

Cleaning Gold: Use an ethanol saturated kimwipe to gently wipe the gold surface thus removing contaminants. Be careful to ensure that no contact is made with the plastic surrounding the gold surface of Spreeta.

Cross-linked Film:

1. Prepare 5 mg/ml bovine serum albumin (BSA) in phosphate buffer saline (PBS), pH 7.4.
2. Prepare 10 mls of 0.4 M EDC and 10 mls 0.1 M NHS, aliquot each into 200 μ l volumes and store at -20 °C.
3. Mix 200 μ l of the EDC preparation with 200 μ l of NHS preparation.
4. Add 0.5 mls of the BSA solution to the EDC/NHS mixture and incubate for 5 min at room temperature.
5. Place 10 μ l the activated BSA solution onto the cleaned gold surface. Spread evenly over the surface. Ensure that the plane of the surface is horizontal and incubate for 15 min at room temperature.
6. Rinse with 20 mM HCl and then in water. Blow dry surface under dry air or nitrogen stream.
7. Add 0.2 g of succinic anhydride to 0.5 mls of DMSO and allow to dissolve (vortexing or sonication will help). Make up to a final volume of 10 mls with 50 mM disodium hydrogen phosphate. Add a few drops of 5M NaOH to adjust pH to around 6.5.
8. Incubate the BSA coated slide in the succinic anhydride preparation overnight at room temperature. Ensure that the gold surface is not in contact with any solid surface. Once coated any physical contact with the gold surface will destroy the film.

Ligand Coupling

Before activating the carboxylated BSA coated surface it is a good idea to test that the surface is properly formed by simply observing the charge preconcentration effect.

Preconcentration Test

1. Dock the sensor with fluidics, initialize, normalize and then equilibrate in water.
2. Dissolve ligand at 20 to 100 µg/ml in 10 mM sodium acetate buffer.
(Note: when preparing the acetate buffer adjust the pH to 1 pH unit below the isoelectric point of the ligand (use 10% acetic acid to adjust pH)). The pH should not be set lower than 4.0 due to protonation of the surface carboxyl groups. Any acidic ligands with pI values less than 4.5 should be immobilized via avidin-biotin linkage (see Application Note 2).
3. Inject 50 µl of the ligand solution at 30 µl /min (ensure that the sensor and all lines are equilibrated in water). Should observe a rapid response due to attraction of the positively charged protein to the negatively charged surface.
4. Rinse with PBS. The high salt concentration in PBS negates the charge interaction and should remove all preconcentrated protein.

Ligand Coupling

1. Activate the surface by injecting a 1:1 mixture of 0.4 M EDC/0.1 M NHS. Allow a surface contact time of 2 - 7 min. Long contact times may convert all the carboxyl groups at the surface to NHS esters and this will destroy the preconcentration effect. It is vital that the majority of carboxyl groups are available to attract the positively charged ligand to the surface. Therefore some trial and error may be required to find the optimum concentration of reagents and appropriate contact time. In some cases EDC/NHS concentrations at 20% of the recommended concentrations proved appropriate.
2. Inject the ligand solution (prepared during the preconcentration test) for 10 min at room temperature. Should observe a response due to both preconcentration and coupling.
3. Rinse with water and then inject 1 M ethanolamine, pH 8.5, for 7 min, to cap residual NHS esters. It is preferable to use PBS, or HBS buffer, rather than water for continuous flow but one must ensure that dispersion within the system is low when using high salt buffer.
4. Check non-specific binding (NSB) using BSA and Ovalbumin at 1 mg/ml in PBS. This also doubles as a blocking step.

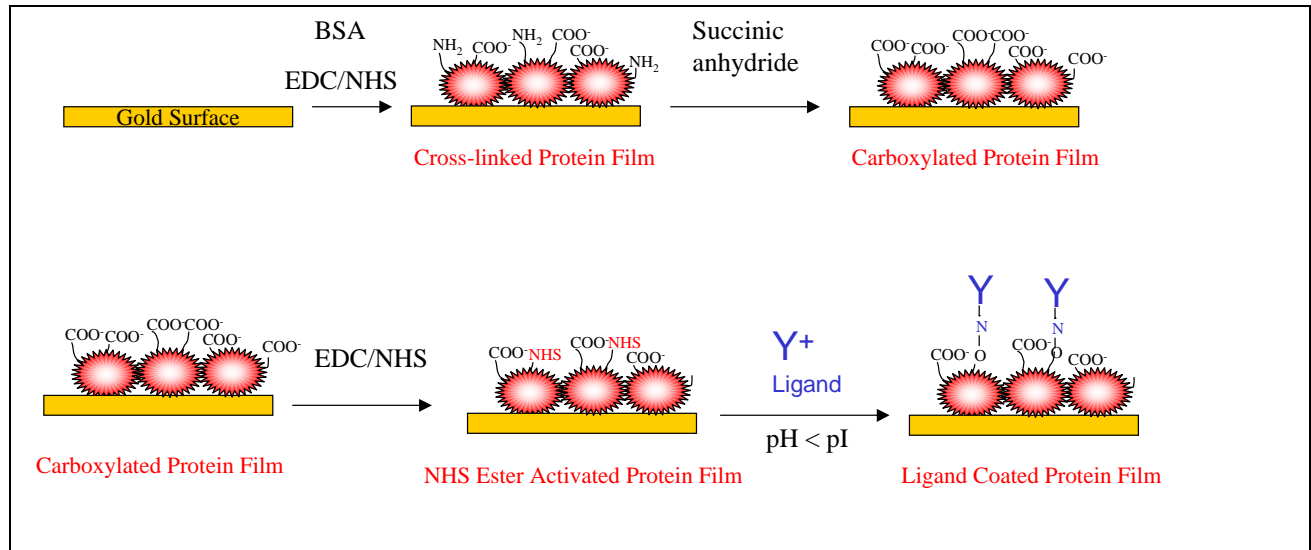


Figure 1. Preparation of cross-linked protein film and immobilization of Ligand

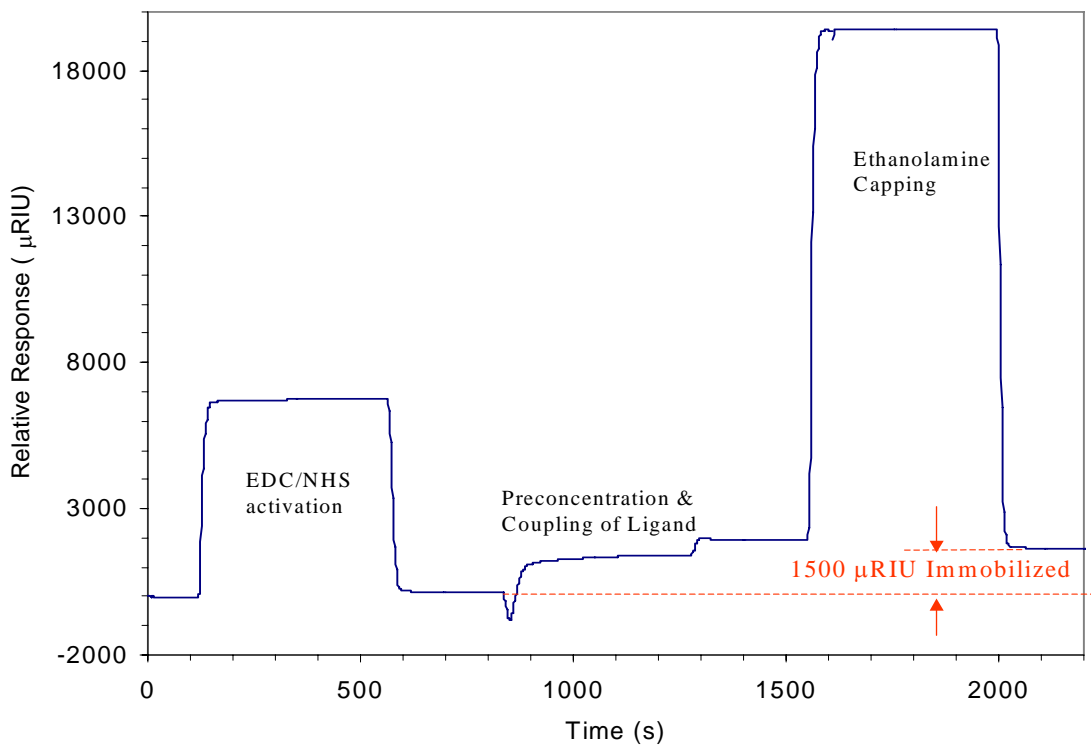


Figure 2. Response curve for immobilization of Goat IgG onto carboxylated protein film by amine coupling. PBS was employed as running buffer.

